


SHINWA EXPRESS	<i>Application</i>	2019
		J u n e
		VOL. 155

Analysis of melamine and related substances using the ULTRON AF-HILIC-CD (HT)

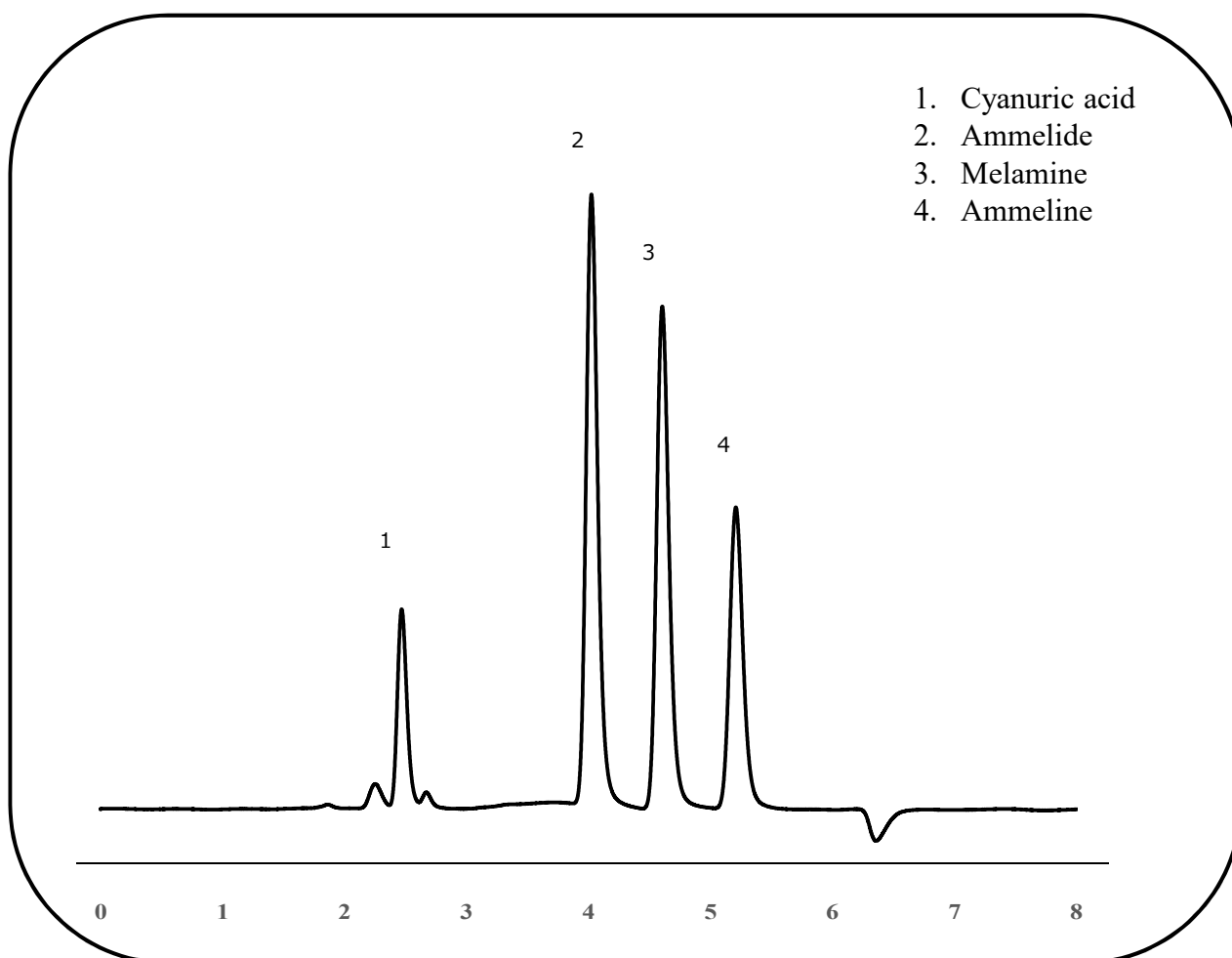

Column: ULTRON AF-HILIC-CD (HT)
Column size: 100 mm L x 2.0 mm I.D.
Mobile phase: 20 mM HCO₂NH₄ (pH 6.8)/CH₃CN = 18/82
Flow rate: 0.15 mL/min

Column temp.: 30°C
Detection: UV-220 nm
Injection vol.: 0.8 µL

●Features of the ULTRON AF-HILIC-CD

1. ULTRON AF-HILIC-CD is a HILIC column in which β -cyclodextrin is chemically bonded to a silica gel via a spacer.
2. The column has high selectivity for molecular structure.
3. The column shows high resolution analysis for structural analogues and structural isomers.
4. The column is suitable for sugar analysis by suppressing schiff base formation and anomer separation.



SHINWA CHEMICAL INDUSTRIES LTD.

50-2 Kagekatsu-cho, Fushimi-ku
Kyoto 612-8307 Japan
TEL: +81-75-621-2360 FAX: +81-75-602-2660
E-mail: info@shinwa-cpc.co.jp
Website: <http://shinwa-cpc.co.jp/en/>

■ Analytical column specifications ■

Product name	Particle size (μm)	Column dimensions Length × I.D. (mm)
ULTRON AF-HILIC-CD (HT)	2	50 × 2.0
		100 × 2.0
		150 × 2.0
ULTRON AF-HILIC-CD	5	100 × 2.0
		150 × 2.0
		100 × 4.6
		150 × 4.6
		250 × 4.6
ULTRON AF-HILIC-DA	5	100 × 2.0
		150 × 2.0
		100 × 4.6
		150 × 4.6
		250 × 4.6

※ Please do not hesitate to contact us for the other dimensions.



Please feel free to contact us with questions related to analyses.
Column screening services are also available.

Please be aware that specifications and prices are subject to change without prior notification.