

SHINWA EXPRESS

Application

2019
August
VOL. 159

Analysis of underivatized amino acids using the ULTRON AF-HILIC-CD (HT-PLSS*)

* PLSS: PEEK-Lined stainless steel column

Column: ULTRON AF-HILIC-CD (HT-PLSS) (2 μ m)

Column size: 150 mm L x 2.0 mm I.D

Mobile phase A: 100 mM HCO₂NH₄ (pH 3.0)/CH₃CN = 10/90Mobile phase B: 100 mM HCO₂NH₄ (pH 3.0)/CH₃CN = 70/30

Flow rate: 0.36 mL/min


Gradient: 20 minutes gradient, 88% - 60% with 10 minutes re-equilibration

Time (min)	%A	%B
0.00	96	4
20.00	50	50
20.01	96	4
30.00	96	4

Column temp.: 40°C

Injection vol.: 1.0 μ LSamples: 5.0 μ g/mL, 5 ppm each

Detection: LC/MS/MS (ESI, Positive, MRM)

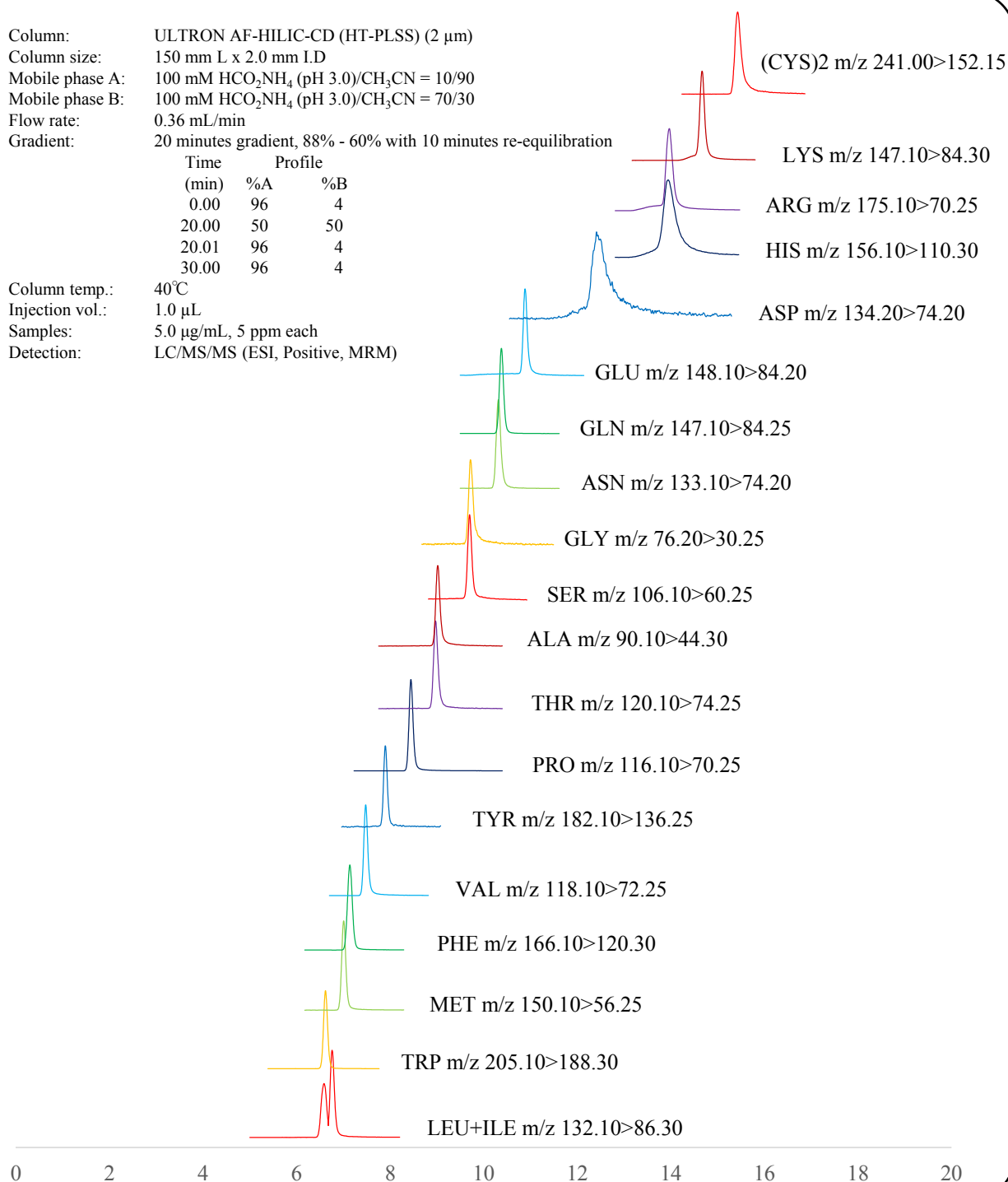

SHINWA CHEMICAL INDUSTRIES LTD.

50-2 Kagekatsu-cho, Fushimi-ku
Kyoto 612-8307 Japan
TEL: +81-75-621-2360 FAX: +81-75-602-2660
E-mail: info@shinwa-cpc.co.jp
Website: <http://shinwa-cpc.co.jp/en/>

● Features of the ULTRON AF-HILIC-CD

1. ULTRON AF-HILIC-CD is a HILIC column in which β -cyclodextrin is chemically bonded to a silica gel via a spacer.
2. The column has high selectivity for molecular structure.
3. The column shows high resolution analysis for structural analogues and structural isomers.
4. The column is suitable for sugar analysis by suppressing schiff base formation and anomer separation.

■ Analytical column specifications ■

ULTRON AF-HILIC-CD (HT)

Product name	Particle size (μm)	Column dimensions Length × I.D. (mm)
ULTRON AF-HILIC-CD (HT)	2	50 × 2.0
		100 × 2.0
		150 × 2.0
		50 × 3.0
		100 × 3.0
		150 × 3.0
ULTRON AF-HILIC-CD (HT-PLSS*)		50 × 2.0
		100 × 2.0
		150 × 2.0
ULTRON AF-HILIC-CD (HT) (Two Guard Cartridges)		
Holder for Guard Cartridge (with Adaptor)	For 5 x 2.0 mm column	

* PLSS: PEEK-Lined stainless steel column

ULTRON AF-HILIC-CD

ULTRON AF-HILIC-CD		
Product name	Particle size (μm)	Column dimensions Length × I.D. (mm)
ULTRON AF-HILIC-CD	5	100 × 2.0
		150 × 2.0
		100 × 4.6
		150 × 4.6
		250 × 4.6
ULTRON AF-HILIC-CD.G		10 × 4.0
ULTRON AF-HILIC-CD (Two Guard Cartridges)		5 × 2.0
		10 × 4.6
Holder for Guard Cartridge (with Adaptor)	For 5 x 2.0 mm column	
	For 10 x 4.6 mm column	

ULTRON AF-HILIC-DA

ULTRON AF-HILIC-DA		
Product name	Particle size (μm)	Column dimensions Length × I.D. (mm)
ULTRON AF-HILIC-DA	5	100 × 2.0
		150 × 2.0
		100 × 4.6
		150 × 4.6
		250 × 4.6
5 × 2.0		
10 × 4.6		
ULTRON AF-HILIC-DA (Two Guard Cartridges)		
Holder for Guard Cartridge (with Adaptor)	For 5 x 2.0 mm column	
	For 10 x 4.6 mm column	

※ Please do not hesitate to contact us for the other dimensions.

Please feel free to contact us with questions related to analyses.
Column screening services are also available.

Please be aware that specifications and prices are subject to change
without prior notification.