


<h1>SHINWA EXPRESS</h1>	<i>Application</i>	VOL. 188
-------------------------	--------------------	----------

Analysis of lower alcohols using Sunpak-A

Sunpak-A is an ethylvinylbenzene-divinylbenzene polymer packing material.

Sunpak-A 80/100

1. Methanol
2. Acetaldehyde
3. Ethanol
4. iso-Propanol
5. n-Propanol
6. sec-Butanol
7. iso-Butanol
8. n-Butanol

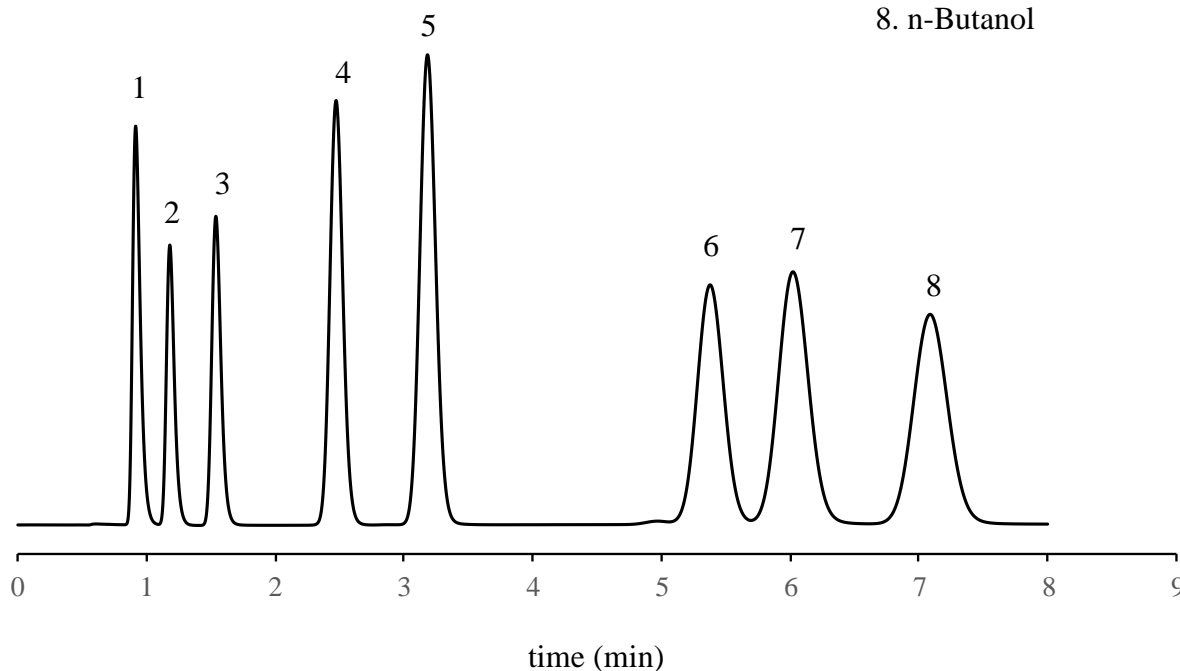

Column: 2.0 m x 3.0 mm I.D. Glass
 Column temp.: 170°C
 Injection port temp.: 220°C

Detector: FID 220°C
 Carrier gas: 50 mL/min N₂
 Sample vol.: 1.0 µL



SHINWA CHEMICAL INDUSTRIES LTD.

50-2 Kagekatsu-cho, Fushimi-ku
 Kyoto 612-8307 Japan
 TEL: +81-75-621-2360 FAX: +81-75-602-2660
 E-mail: info@shinwa-cpc.co.jp
 Website: <https://shinwa-cpc.co.jp/en/>