


SHINWA EXPRESS


Application

VOL. 191


Analysis of Halogenated acetic acid using the ULTRON PS-80H

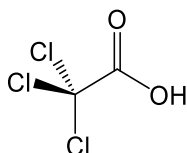


1. Trichloroacetic acid

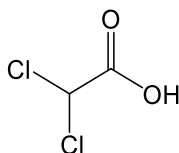


3. Dichloroacetic acid

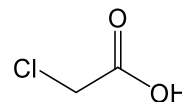


5. Chloroacetic acid

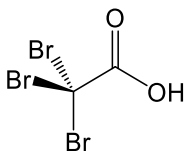


2. Tribromoacetic acid

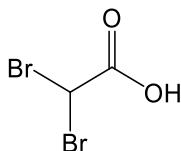

4. Dibromoacetic acid

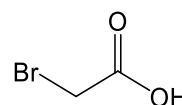

6. Bromoacetic acid

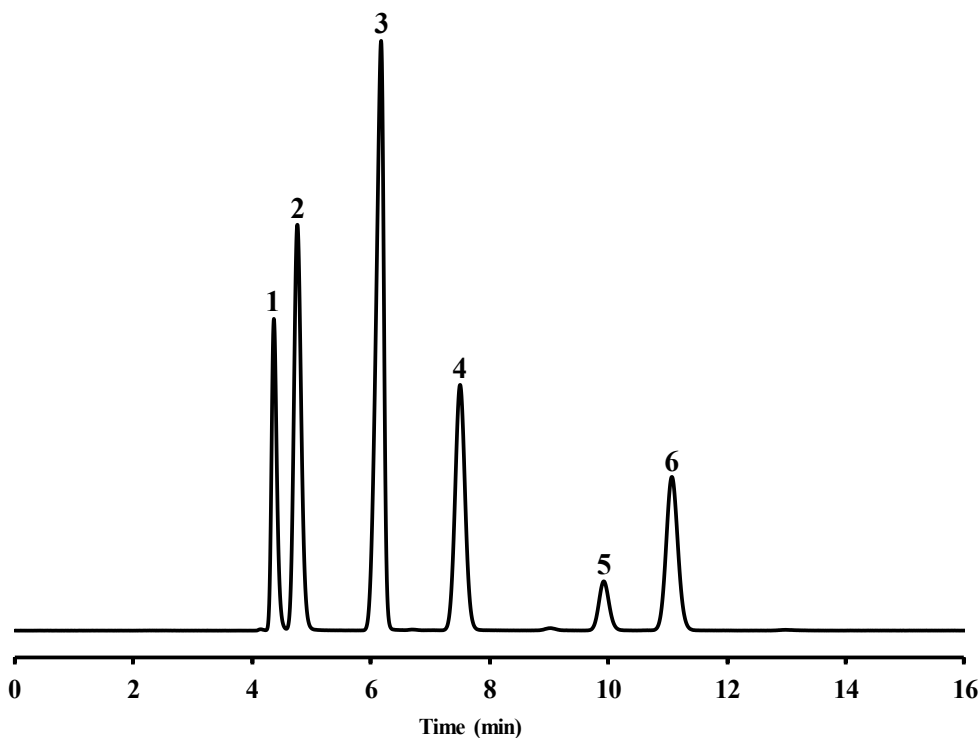

Column: ULTRON PS-80H (10 μ m)
 Column size: 300 mmL x 8.0 mm I.D.
 Mobile phase : HClO_4 aq (pH 2.1)
 Flow rate: 1.0 mL/min

Column temp.: 60°C
 Detection: UV-210 nm
 Injection vol.: 10 μ L



SHINWA CHEMICAL INDUSTRIES LTD.

50-2 Kagekatsu-cho, Fushimi-ku
 Kyoto 612-8307 Japan
 TEL: +81-75-621-2360 FAX: +81-75-602-2660
 E-mail: info@shinwa-cpc.co.jp
 Website: <https://shinwa-cpc.co.jp/en/>

● Features of the ULTRON PS-80H

1. The ULTRON PS-80H columns are HPLC columns packed with a sulfonic acid type polystyrene cation exchange resin designed for organic acid analyses.
2. The columns utilize electrostatic repulsion interactions.
3. The columns display excellent retention reproducibility.
4. The columns are durable due to the use of a mechanically strong polymer gel.

■ Analytical column specifications ■

ULTRON PS-80H / PS-80N / PS-80C / PS-80P

Product name	Particle size (μm)	Column dimensions Length x I.D. (mm)
ULTRON PS-80H	10	300×8.0
		250×2.0
ULTRON PS-80H.G		50×8.0
ULTRON PS-80N		300×8.0
ULTRON PS-80N.L		500×8.0
ULTRON PS-80N.G		50×8.0
ULTRON PS-80C		300×8.0
ULTRON PS-80C.G		50×8.0
ULTRON PS-80P		300×8.0
ULTRON PS-80P.G		50×8.0



※ Please do not hesitate to contact us for the other dimensions.
Please feel free to contact us with questions related to analyses.
Column screening services are also available.

Please be aware that specifications and prices are subject to change without prior notification.