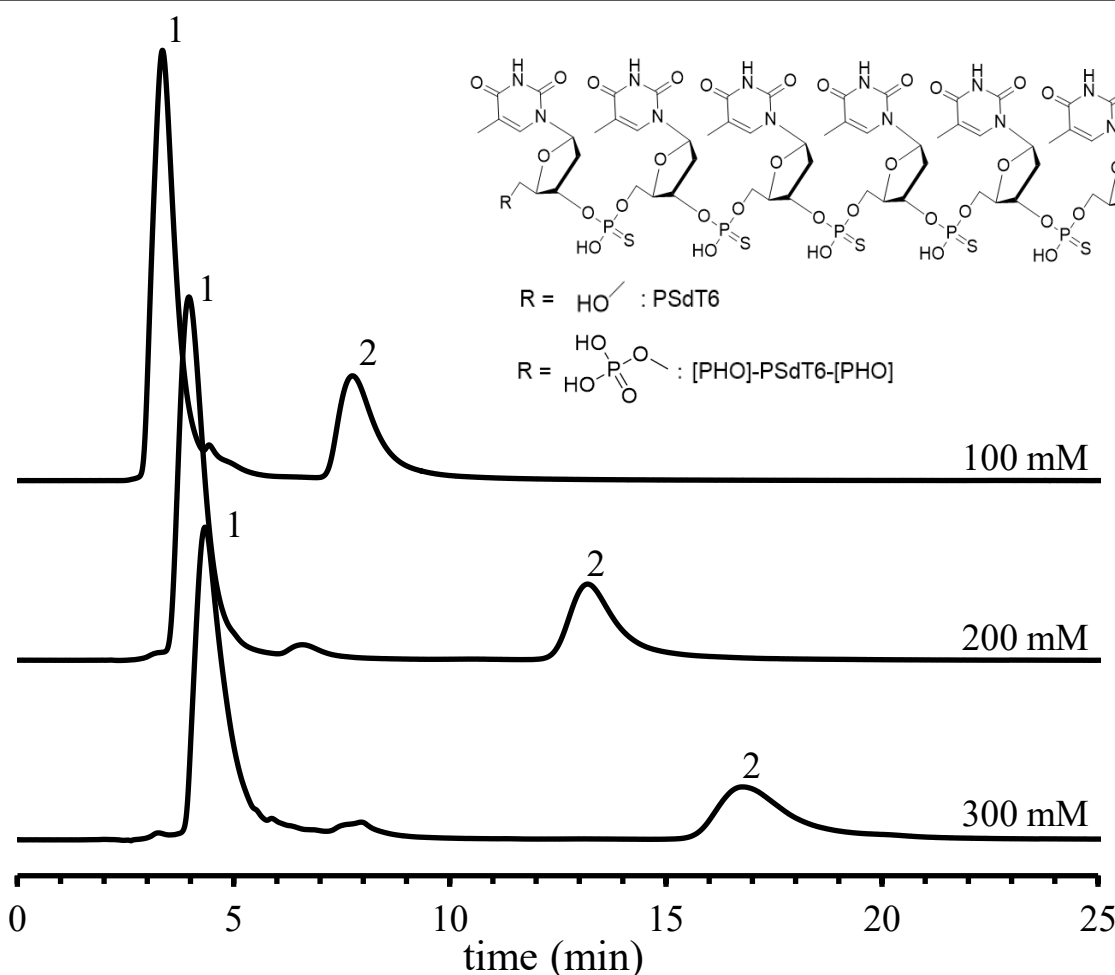


SHINWA EXPRESS

Application

VOL. 256

Analysis of PSdT6 and [PHO]-PSdT6-[PHO] using the ULTRON AF-HILIC-CD (Part 1) “Effect of salt concentration in mobile phase”

Column: ULTRON AF-HILIC-CD (2 μm)

Column size: 150 mmL x 2.0 mm I.D.

Mobile phase: 100 mM, 200 mM, 300 mM $\text{CH}_3\text{CO}_2\text{NH}_4/\text{CH}_3\text{CN} = 35/65$

Flow rate: 0.15 mL/min

Column temp.: 40°C

Detection: UV-260 nm

Sample: 1. 5'-T*T*T*T*T*T-T-3'

2. 5'-[PHO]-T*T*T*T*T*T-T-[PHO]-3'

Diluent: $\text{CH}_3\text{CN}/\text{TE buffer}^* = 70/30$

* TE buffer = 10 mM Tris-HCl buffer pH 8.0 +

1 mM EDTA buffer pH 8.0

Injection vol.: 0.5 μL

● Features of the ULTRON AF-HILIC-CD

1. ULTRON AF-HILIC-CD is a HILIC column in which β -cyclodextrin is chemically bonded to a silica gel via a spacer.
2. The column has high selectivity for molecular structure.
3. The column shows high resolution analysis for structural analogues and structural isomers.
4. The column is suitable for sugar analysis by suppressing schiff base formation and anomer separation.



SHINWA CHEMICAL INDUSTRIES LTD.

50-2 Kagekatsu-cho, Fushimi-ku
Kyoto 612-8307 Japan

TEL: +81-75-621-2360 FAX: +81-75-602-2660

E-mail: info@shinwa-cpc.co.jp

Website: <https://shinwa-cpc.co.jp/en/>

■ Analytical column specifications ■

ULTRON AF-HILIC-CD

Product name	Particle size (μm)	Column dimensions Length x I.D. (mm)
ULTRON AF-HILIC-CD	2	50×2.0
		50×3.0
		100×2.0
		100×3.0
		150×2.0
		150×3.0
ULTRON AF-HILIC-CD (PLSS)※ ※ PLSS: PEEK-Lined Stainless steel		50×2.0
100×2.0		
150×2.0		
ULTRON AF-HILIC-CD (Two Guard Cartridges)		5×2.0
ULTRON AF-HILIC-CD	5	100×2.0
		100×4.6
		150×2.0
		150×4.6
		250×2.0
		250×4.6
ULTRON AF-HILIC-CD.G		10×4.0
ULTRON AF-HILIC-CD (Two Guard Cartridges)		5×2.0
		10×4.6
Holder for Guard Cartridge (with Coupler)		-
	For 10×4.6 mm column	

ULTRON AF-HILIC-DA

Product name	Particle size (μm)	Column dimensions Length x I.D. (mm)
ULTRON AF-HILIC-DA	5	100×2.0
		100×4.6
		150×2.0
		150×4.6
		250×4.6
		5×2.0
ULTRON AF-HILIC-DA (Two Guard Cartridges)		10×4.6
Holder for Guard Cartridge (with Coupler)	-	For 5×2.0 mm column
		For 10×4.6 mm column



※ Please do not hesitate to contact us for the other dimensions.
Please feel free to contact us with questions related to analyses.
Column screening services are also available.

Please be aware that specifications and prices are subject to change without prior notification.