

# WCOT Capillary Column

## For Gas Chromatography

### INSTRUCTION MANUAL

#### 1. Caution

Please read these instructions carefully before using this product.

#### 2. Instruction

- ULBON WCOT capillary columns are columns in which liquid phases are coated on internal walls. Use this column only for gas chromatography.

#### 3. Notes

- As the retention of stationary phases on open tubular columns (WCOT, SCOT) is weaker than that of packed columns, these columns are susceptible to rapid temperature changes with the result that rapid changes in temperature can lead to the deterioration of these columns.

Consequently, when raising the temperature of this column to the intended analytical temperature and when using temperature programming, please ensure that the rate of increase is 6°C/min or less.

The maximum operating temperature for this column is 160°C.



50-2 Kagekatsu-cho, Fushimi-ku, Kyoto 612-8307 JAPAN

TEL 075-621-2360 URL <https://shinwa-cpc.co.jp>